

# medilines

# Albert's Colour

For Diphtheria Bacilli

## REAGENTS & PREPARATIONS

ALBERT'S SOLUTION -1 : Ready to use.  
ALBERT'S SOLUTION -2 : Ready to use.

## PREPARATION & FIXATION OF SMEAR

Fix the dry smear by heating on slide.

## STAINING PROCEDURE

- Cover the fixed slide with **ALBERT SOLUTION - 1** for 2 minutes.
- Rinse with Water and blot dry.
- Counter stain with **ALBERT SOLUTION - 2** for 1 minute.
- Rinse with Water, Blot dry & Examine under the microscope.

## RESULT

Granules of Diphtheria Bacilli	Black
Cytoplasm	Light green