

medilines

Fungus Colour

For Fungus staining

REAGENTS & PREPARATIONS

medilines Fungus Colour -Ready to use

STAINING PROCEDURE

1. Place a drop of 95% alcohol on the slide.
2. Gently tease out a fragment of the culture in the alcohol with needle or straight wire. When it is satisfactorily spread, let most of the alcohol evaporates.
3. Add a drop of medilines Fungus Colour.
4. Apply a cover glass, avoiding bubbles, and exert gentle pressure if the fungus fragments do not lie flat.
5. Remove any excess stain round the cover glass with the edge of a blotting paper.
6. Let the stain penetrate. (It may be satisfactory within a few minutes but differentiation may go on improving for up to 24 hours.)
7. For permanent preparations, seal the edge of cover glass with nail varnish or cellulose lacquer.
8. View under a microscope.

medilines

medilines@medilines.com