

medilines

H & E COLOUR

HAEMATOXYLIN AND EOSIN

REAGENTS & PREPARATIONS

medilines Haematoxylin:	Ready to use
medilines Decolourizer:	Ready to use
medilines Eosin:	Ready to use

STAINING PROCEDURES

1. Place **medilines** Hematoxylin on the specimen for 4-5 minutes
2. Rinse in distilled water.
3. Place in **medilines** Decolouriser for a few seconds
4. Rinse in distilled water.
5. Place **medilines** Eosin on the specimen for counter staining
6. Wash in tap water

Note

- The above Staining Procedure is a routine procedure. For specific applications the procedure can be modified by yourself
- In certain procedures, add 2-4 ml Acetic acid per 100ml Haematoxylin. It is optional but gives sharper nuclear staining. this must be added just before use.
- Blueing Solutions and Mounting medium are also available.